

Initial Release July, 2024
Revised January, 2025

TECHNICAL PRODUCT BULLETIN TB-2024-07-01 Rev B

Subject: LD Series Accessory Changes

25LD Tipping Stand


As previously detailed in Technical Bulletin TB-2024-03-01 there have been changes to the LD Series product line with the 25LDB vessel becoming obsolete and replaced by the 25LD. Due to this change the tipping stand, item number D025-8C00 is not compatible with the new 25LD.

- If you require a tipping stand for the older 25LDB with a 15.6" (396mm) outer diameter then please order the original tipping stand, item number D025-8C00.
- If you require a tipping stand for the new 25LD with a 14.8" (376mm) outer diameter then please order the new version tipping stand, item number D025-8D01.

Liquid Withdrawal Device

The liquid withdrawal device for the 25LD, 35LD and 50LD has been changed so it is compatible with both the old design LDB vessels and the new design LD vessels. Previously the liquid withdrawal tube of the liquid withdrawal device had a single score approximately 6 in. (152 mm) from the end. When using the withdrawal device with the 25 and 35 liter vessels the scored section was broken off to make the tube shorter.

Now the liquid withdrawal tube of the liquid withdrawal device has three score marks as shown in the drawing below.

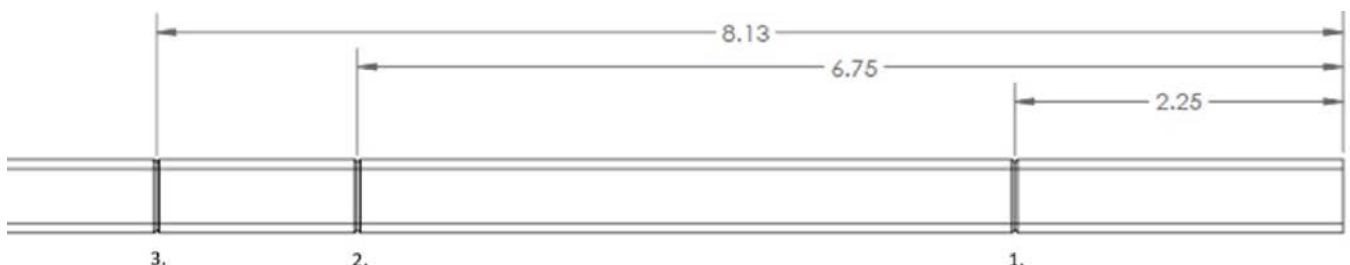

 1-855-750-8191

 www.icbiomedical.com

 1031 Cass White Road, NW
Cartersville, Georgia 30121



- The liquid withdrawal tube without breaking off any of the sections is for use with the older 50LDB vessel.
- Breaking off the tube at point 1. the first score from the bottom of the tube is for use with the newer 50LD vessel.
- Breaking off the tube at point 2. the second score from the bottom of the tube is for use with the 35LD, 35LDB and the older 25LDB vessels.
- Breaking off the tube at point 3. the third score from the bottom of the tube is for use with the newer 25LD vessels.

For customers who have already purchased older versions of the liquid withdrawal device with the single score mark withdrawal tube and wish to use it with the newer 25LD. You have the choice to either cut the existing withdrawal tube by 1.38" (35mm) or you can purchase a new withdrawal tube (item number D050-8C22) that has the 3 score lines.

 1-855-750-8191

 www.icbiomedical.com

 1031 Cass White Road, NW
Cartersville, Georgia 30121